

Software 4.0 upgrade guide



 **SERSTECH**
Chemical intelligence solutions ■

Software upgrade guide – to version 4.0

Make a backup first!

-  Your old measurements and local databases will not automatically follow the upgrade. If you want to keep your history and local databases, please remember to make a backup before you proceed with the upgrade.


How to make a backup

- Connect the instrument to a PC using a USB cable
- Use the quick menu to activate USB.




How to make a backup

➤ Activate USB DATA

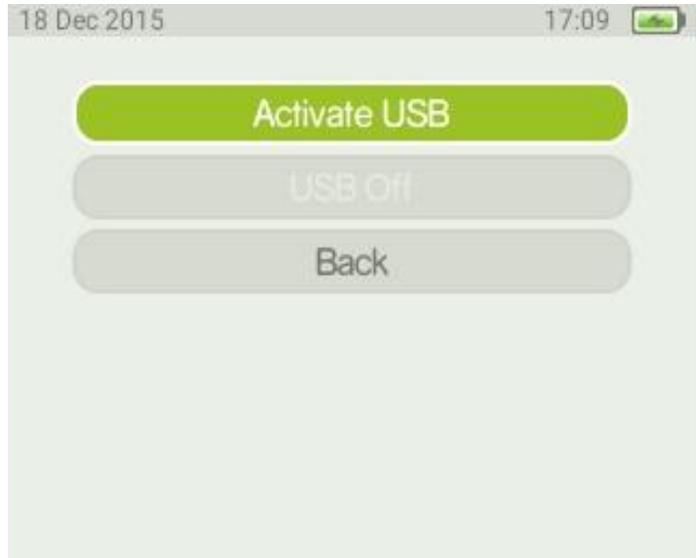


➤ A unit called i100_data will appear on your computer

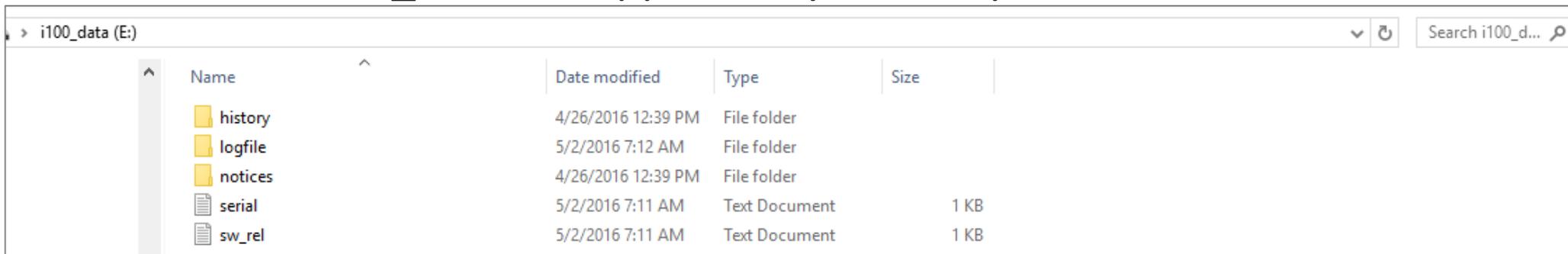

How to make a backup

- Open i100_data and copy the history folder. Open i100_conf and copy your local databases. Save it to a location on your PC or other location outside the Indicator.
- Please note, once you have made the upgrade to v.4.0 you cannot directly put the files back on the Indicator, you will need to use ChemDash. Contact Serstech Support, support@serstech.com for more information.
- Select USB Off on the Indicator.

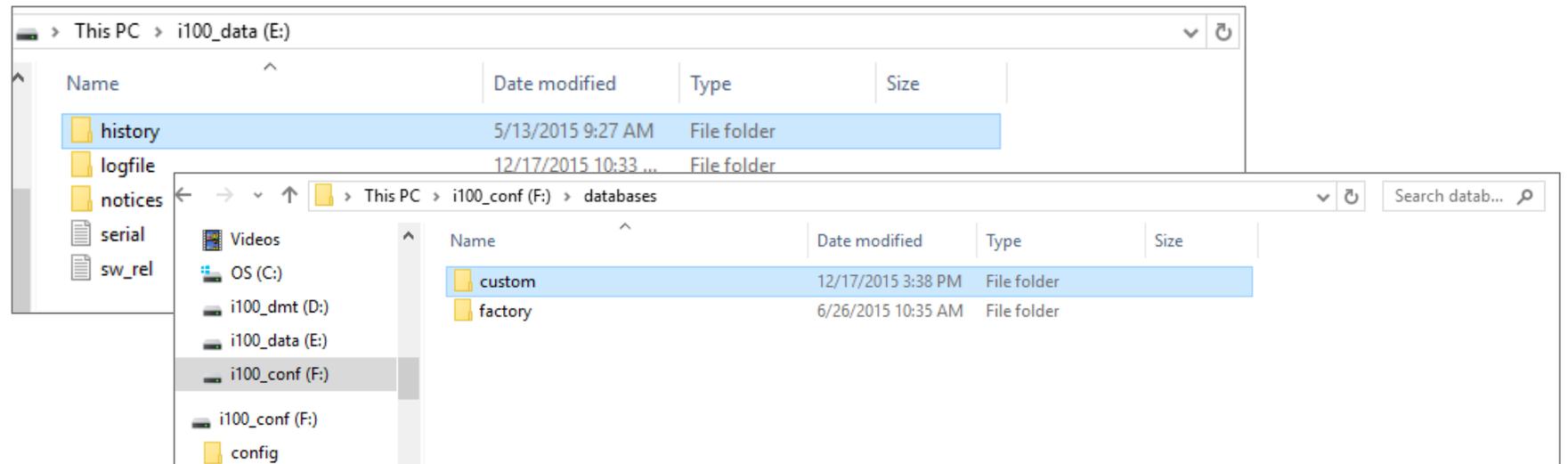

How to upgrade to 4.0

- First, make sure the battery of your instrument is fully loaded.
- **Please check your current Software version first!** Make sure your instrument has the latest release 3.3.6 or later before upgrading to release 4.0. If the indicator has the wrong software version, the upgrade will fail.
- If you are unsure about your current version, go to “Settings” -> “About” in your Indicator.
- Download the file i100update 4.0 or if you have an earlier version than 3.3.6, make a first upgrade to 3.3.6.
- Save the file on your computer, not on an USB stick, that can sometimes create conflicts.

Name	Date modified	Type	Size
 i100update-REL-4.0.1-r2827-20160429-13...	4/29/2016 1:36 PM	GZ File	42,665 KB


Upgrade using "Drag and Drop"

- ⇨ Log in to your Indicator and activate its USB mode:
 - ⇨ Settings->Administration->USB->Activate USB
 - ⇨ Or use "Quick Menu" ->USB-> Activate USB
 - ⇨ Note! You need to have the read and write privileges to do the updates. This can be set by the admin user.
- ⇨ Connect the Indicator to your computer.
- ⇨ Look for the i100_conf in your file browser



Upgrade using "Drag and Drop"

- We strongly recommend that you read the Release Notes before upgrading.
- Drag and drop the file to the folder "**updates**" in your instrument (i100_Conf)

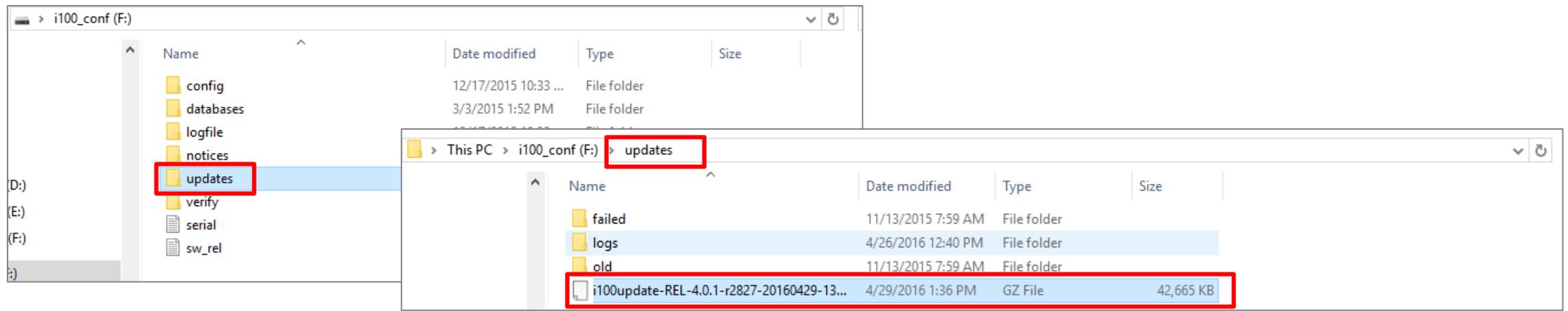

- Deactivate USB, unplug from your computer and restart the instrument. Restarting can take time (up to 15 minutes). Please be patient.
- To see which Software version you have, turn on your instrument, go to Setting -> About.

Software upgrade guide – to v4.0.

You should now have upgraded successfully!

