

Get started – connect to ChemDash using the network connection kit

Wireless connection - Wifi

■ **1. Please upgrade your Serstech 100 Indicator to software 4.x, otherwise ChemDash cannot be connected.** If your Indicator runs an older version please upgrade the software first by following the instructions in the Software Upgrade Guide on serstech.com.

■ The Network connection kit content



USB Y-cable for dongle connection

To Indicator

To WiFi

To power supply



Ethernet dongle



WiFi dongle


■ 2. Connect the Indicator to your PC with the delivered USB cable



USB cable



■ 3. Login to the Indicator

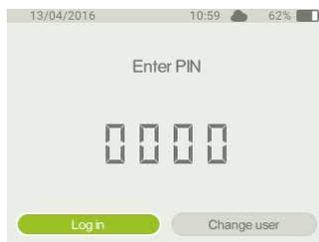

■ 4. Choose USB in the quick menu on the Indicator Keypad and activate USB



Get started – connect to ChemDash using the connection kit

Wireless connection – WiFi


5. Go to your PC



5.1 Run the wifi settings tool from the Indicator

5.2 Choose the wireless network you want to use from the list and enter the password. Click update.

6. Connect USB Y-cable as shown

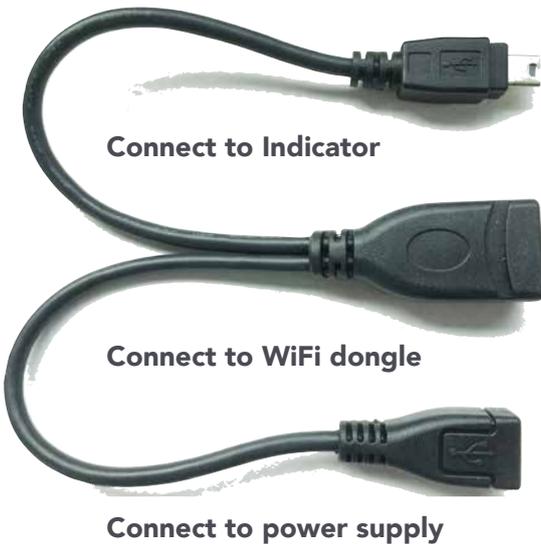

Indicator



WiFi dongle



Micro USB cable + 1A USB power adapter (not delivered in the case)

7. Go to www.chemdash.com and login with the login default you have received by email from Serstech.

Find your device by it's serial number (shown on the backside of the instrument) in the Device Management with a green online status.

Device	Device Tag	Status	Last Seen	Model	Software Version	Calibration D
000000000060		Online	2016-05-23 14:05	Serstech 100 Indicator	4.0.2	2016-05-23 01:...
000000000089		Offline	2016-05-23 12:43	Serstech 100 Indicator	4.0.2	2016-04-25 12:33

